

Template upload form

Name of Template	ControlVaccumBox BVSVaccum
Name of designer	Martin Zogg
Purpose of the Template	<p>This is a template to regulate the vacuum pressure in the vacuum box.</p> <p>Two thresholds can be specified:</p> <ul style="list-style-type: none"> - the end pressure that shall be reached - a threshold pressure (after most of the liquid is transferred). <p>The system will return the information if only the threshold pressure, both threshold and target pressure or none of the pressures has been reached.</p> <p>If the one of the pressures is not reached, a user output is shown. The user then has the possibility to press the button Retry for a second run, Ignore to go forwards or Cancel to abort.</p> <p>Please note that the pressure must be entered as delta pressure, so it is not an absolute value but a difference to the ambient pressure.</p>
Creation date	February 19, 2008
Tested on	
Vector software version	4.2.0.6263
Service Pack of Vector software	-
Operating system	Windows XP professional Version 2002
Service Pack(s) of OS system	Service Pack 2
Instrument	ML StarLet SN 1470
Instrument Configuration	BVS Pump, Carrier 190032/00
Firmware version(s)	-
Parameters to pass	<ul style="list-style-type: none"> - i_pumpId - i_comPort - i_simulatorMode - i_vacDeltaPressure - i_vacThreshold - i_vacTime - i_vacValveOpen
Search strings	
Functions	-
How to install	<p>Save the template files in the subfolder "Templates". Using templates for the first time, the folder "Templates" has to be created. The correct spelling of this folder name is important:</p> <p>C:/Program files/Hamilton/Library/Templates</p>
How to uninstall	<p>Delete Template files in the folder:</p> <p>(C:/Program files/Hamilton/Library/Templates)</p>

Template upload form

Name of Template	ControlVaccumBoxBVSAbsort
Name of designer	Martin Zogg
Purpose of the Template	This template it used to de-Init the com interface
Creation date	19. February 2008
Tested on	
Vector software version	4.2.0.6263
Service Pack of Vector software	-
Operating system	Windows XP professional Version 2002
Service Pack(s) of OS system	Service Pack 2
Instrument	ML StarLet SN 1470
Instrument Configuration	BVS Pump, Carrier 190032/00
Firmware version(s)	-
Parameters to pass	- i_pumpId - i_comPort
Search strings	
Functions	-
How to install	Save the template files in the subfolder "Templates". Using templates for the first time, the folder "Templates" has to be created. The correct spelling of this folder name is important: C:/Program files/Hamilton/Library/Templates
How to uninstall	Delete Template files in the folder: (C:/Program files/Hamilton/Library/Templates)